

Technical Procedure for Ardrox

1.0 Purpose – This procedure is a non-porous development procedure.

2.0 Scope – This procedure is used in the processing of non-porous evidence that may contain impressions that require developing/enhancing.

2.1 One of the most effective ways to recover latent prints from items of evidence is to use a fluorescent dye followed by a laser or alternate light source examination. Ardrox is one of the most effective laser dyes for use on various non-porous surfaces. This dye is normally for use on non-porous surfaces; however, it may, under certain conditions, be used on porous or semi-porous surfaces. Ardrox is extremely efficient; it is highly fluorescent and can be used with various alternate light sources.

3.0 Definitions

- **Alternate light source:** Any of the multiple forensic light sources readily available in the Digital/Latent Evidence Section including, but not limited to, the CrimeScope, Mini-CrimeScope, TracER Laser, and Ultra-Lite ALS.
- **CE:** Cyanoacrylate ester also known as super glue.

4.0 Equipment, Materials and Reagents

4.1 Equipment and Materials

- Orange filter (goggles and/or camera lens filter)
- Alternate light source
- Fume hood
- Gloves
- Face shield and/or safety goggles
- Plastic applicator bottles or tray for submersion

4.2 Reagents

- Ardrox concentrate solution
- Methanol

5.0 Procedure

5.1 Mixing Procedure

5.1.1 Place six (6) mL of Ardrox concentrate solution in one hundred (100) mL of methanol. Stir until solution is completely incorporated into the methanol. Solution is ready for use.

5.2 Application Procedure

5.2.1 Forensic Scientists shall produce a self-made test print to be processed concurrently with items of evidence. (See Section Technical Procedure for Ensuring Quality Control.)

5.2.2 First apply cyanoacrylate ester to the item of evidence. Ardrox adheres to the chlorosis that occurs after processing with CE. Utilizing the fume hood and gloves, spray, immerse, or brush item of evidence with the methanol/Ardrox solution. Allow to air dry for at least one minute.

5.2.3 Rinse item with running tap water to remove any excess Ardrox solution. Allow to air dry (this is especially important for semi-porous items).

5.2.4 When completely dry, view the item using the CrimeScope or Mini-CrimeScope (300 – 400 nm). Wear goggles to view any fluorescence on the item. Latent prints will fluoresce bright yellow.

5.2.5 Any latent prints shall be preserved using photography. Camera shall be equipped with an orange filter for print visualization.

5.3 Standards and Controls – N/A

5.4 Calibration – N/A

5.5 Sampling – N/A

5.6 Calculations – N/A

5.7 Uncertainty of Measurement - N/A

6.0 Limitations – Ardrox is designed to be used for the processing of non-porous evidence. Some porous/semi-porous items may be processed using this technique.

6.1 The cyanoacrylate fuming process is vital to the success of Ardrox processing.

6.2 Ardrox concentrate solution has an indefinite shelf life. Once prepared, it has a shelf life of six (6) months.

6.3 All prepared solution shall be stored in dark, shatter-proof bottles. Working solution may be stored in clear spray bottles.

6.4 Ardrox concentrate solution shall be stored in its original shipping container.

6.5 Ardrox may be used in conjunction with other fluorescent dyes.

6.6 Inherent luminescence of the item to be processed must be conducted before treatment with Ardrox in order to assess if the item fluoresces independently. This will help to determine if Ardrox is the appropriate dye for latent print visualization.

7.0 Safety – The chemical concentrate solution is flammable and must be handled with extreme care. It may cause some irritation when in contact with the skin or eyes and may be harmful if inhaled or ingested. The methanol used in the solution preparation is toxic and must be handled with care. Protective gloves, eye goggles and aprons shall be worn at all times.

8.0 References

Kent, T., ed. *Manual of Fingerprint Development Techniques: A Guide to the Selection and Use of Processing for the Development of Latent Fingerprints*. Police Scientific Development Branch, London (July 1992).

Lee, H.C. "Methods of Latent Print Development." *Proceedings of the International Forensic Symposium on Latent Prints*. (July 1987): 15 – 24.

Lennard, C.J. and P.A. Margot. "Sequencing of Reagents for the Improved Visualization of Latent Fingerprints." *Proceedings of the International Forensic Symposium on Latent Prints*. (July 1987): 141-142.

Manual of Fingerprint Development Techniques. (January 1986): 2-8.

Manual of Fingerprint Development Techniques: A Guide to the Selection and Use of Processes for the Development of Latent Fingerprints. Scientific Research and Development Branch, London (1986).

Olenik, J. "Ardrex: An Alternate Solvent System." *Journal of Forensic Identification*. Vol. 42, 6: 513-516 (1992).

Trozzi, T.A., R.L. Schwartz and M.L. Hollars. *Processing Guide for Developing Latent Prints*. (2000): 1-64.

US Department of Justice. *Chemical Formulas and Processing Guide for Developing Latent Prints*. FBI Laboratory Division, Latent Fingerprint Section (1994).

Wilkinson, D.A., and A.H. Misner. "A Comparison of Thenoyl Europium Chelate with Ardrex and Rhodamine 6G for the Fluorescent Detection of Cyanoacrylate Prints." *Journal of Forensic Identification*. Vol. 44, 4: 387-406 (1994).

9.0 Records - N/A

10.0 Attachments - N/A

Revision History		
Effective Date	Version Number	Reason
09/17/2012	1	Original Document
10/31/2013	2	Added issuing authority to header