

## OVEN CALIBRATION LOG

Date \_\_\_\_\_

## Oven Setting

## Thermometer Reading

Initials

[illegible]

## Trace Evidence

## **Oven**

### **Calibration Procedures**

Trace Evidence ovens will be calibrated monthly. The results of the calibration should be recorded in the oven calibration log.

**Procedure:**

1. Record date.
2. Read oven setting and record in degrees celsius.
3. Read thermometer temperature and record in degrees celsius.
4. Initials should be recorded by chemist conducting calibration.
5. Calibration of oven setting and thermometer should be within +/- 5 degrees celsius.
6. If calibration fails then oven should be recalibrated. (See Note)

Note: For Fisher Scientific Isotemp Ovens see the  
Fisher Scientific Isotemp Instruction Manual

For Blue M Laboratory Oven see the  
Blue M Operation and Maintenance Manual